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
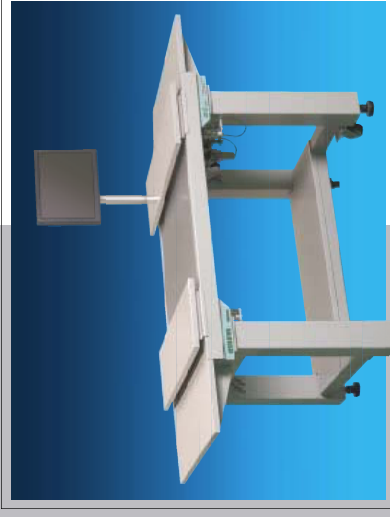
### DESK TYPE AIR SUSPENSION VIBRATION ISOLATOR

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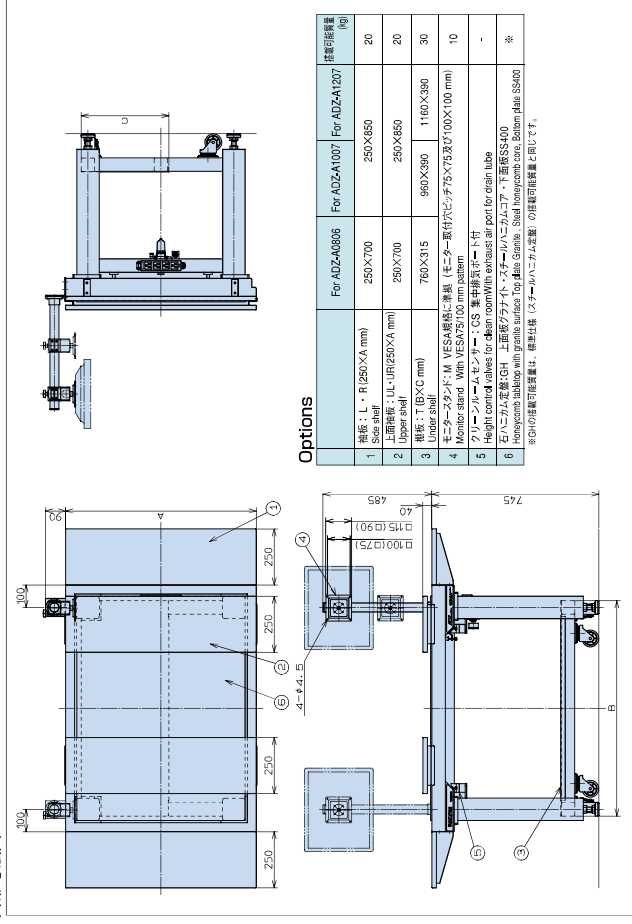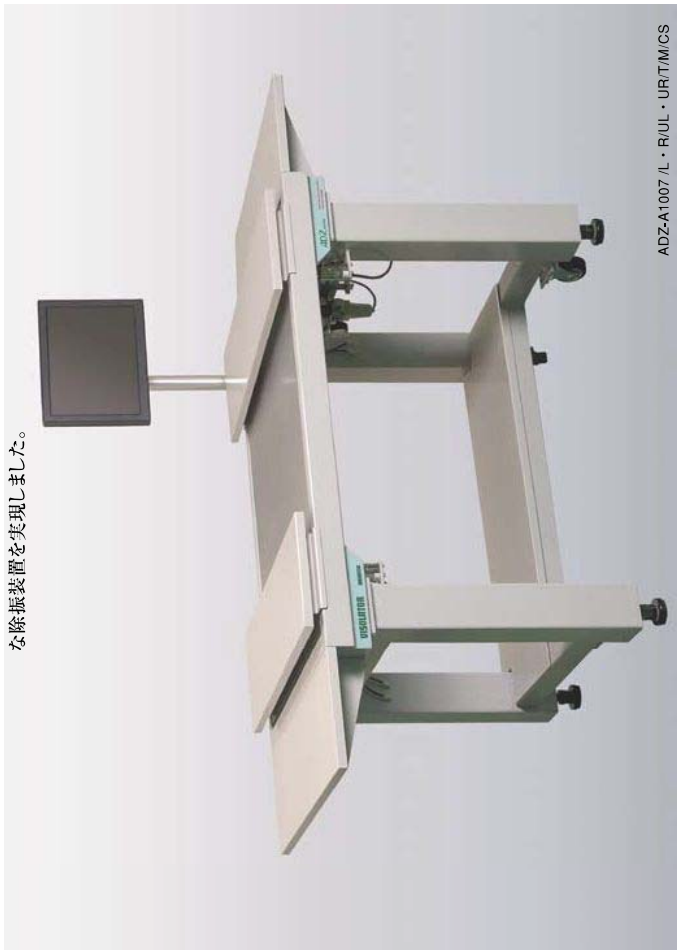
ADZ-A1007/L・R/UL・UR/T/M/CS

The ADZ-A series is a highly upgraded version of the well-selling ADZ series. While its design concept remains the same as that of the ADZ series, it incorporates many new solutions demanded by the customer. Vibration isolation performance is further enhanced, working comfort is greatly improved, and a model with a larger mounting plate has been added to the new series. Also, a wealth of options is available for better operability of the equipment installed on the isolator. This high cost performance vibration isolator can accommodate a wide range of equipment and can be used almost anywhere.

### ■ Has excellent vibration isolation performance in both the vertical and horizontal directions.

The use of a thin diaphragm type air spring, which has a long history and a proven sales track record, together with a new damping rubber, realize enhanced vibration isolation performance in both the vertical and horizontal directions. In addition, cost performance is excellent.

### ■ Design based on the customer's viewpoint

The color and refined exterior, which gives the impression of cleanliness, make this isolator suitable for an office environment as well. The user-friendly design is aimed at more comfortable workability. Three models are available in this series to enable it to cope with a wide range of equipment. Multiple power outlets (four 2-pole sockets + ground terminal) and a urethane armrest are provided as standard fitment.

### ■ Automatic leveling mechanism

The automatic leveling mechanism immediately restores the isolator to a horizontal position after the equipment installed on it is changed over, or the mounting plate becomes inclined due to a shift in the center of gravity. Also, a surge tank and orifice are provided to rapidly dampen shakiness of the mounting plate.

### ■ Use of a steel honeycomb optical tabletop as the mounting plate

This isolator uses a highly rigid, lightweight steel honeycomb optical tabletop which has high dampening performance. The top surface plate, made of SUS430 ferromagnetic stainless steel, enabling a magnet stand to be used. (We can provide tap holes on the top surface, drill holes, and so on, to suit your requirements. Please ask us for details.)

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### ■ A wealth of options

● **Side shelves**：L and R  
These shelves can provide additional space for working or for installing peripheral equipment. They can be installed on the left and right sides of the main unit. Width is 250 mm.

● **Upper shelves**：UL and UR  
These shelves are installed straddling the mounting plate. Because they do not directly touch the mounting plate, they do not impair the vibration isolation performance. These shelves can be slid left and right, and used in any desired position. Width is 250 mm.

● **Under shelf**：T  
The under-shelf can provide storage space beneath the frame.

● **Monitor stand**：M  
The monitor fixing part conforms to the VESA standard (screw pitch: 75×75 mm and 100×100 mm). It enables the height and angle of the monitor to be adjusted. It can be installed at rear right or rear left of the isolator.

● **Height control valve for clean room**：CS  
A drain tube can be connected to this valve to enable the air exhausted from the air spring to be discharged to a suitable location. This enables you to use the isolator in a clean room with peace of mind. (In this case, the standard height control valve is changed to one that has clean room specifications.)

● **Honeycomb tabletop with granite surface**：GH  
Honeycomb structure tabletop with highly corrosion-resistant granite top surface plate

■ **Applications**  
Optical microscopes, laser microscopes, scanning probe microscopes, electrophysiology experiments, microtransmission, interferometers, manual probers, high magnification CCD, surface roughness gauges, others

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