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
## HIGH QUALITY OF VIBRATION ISOLATORS

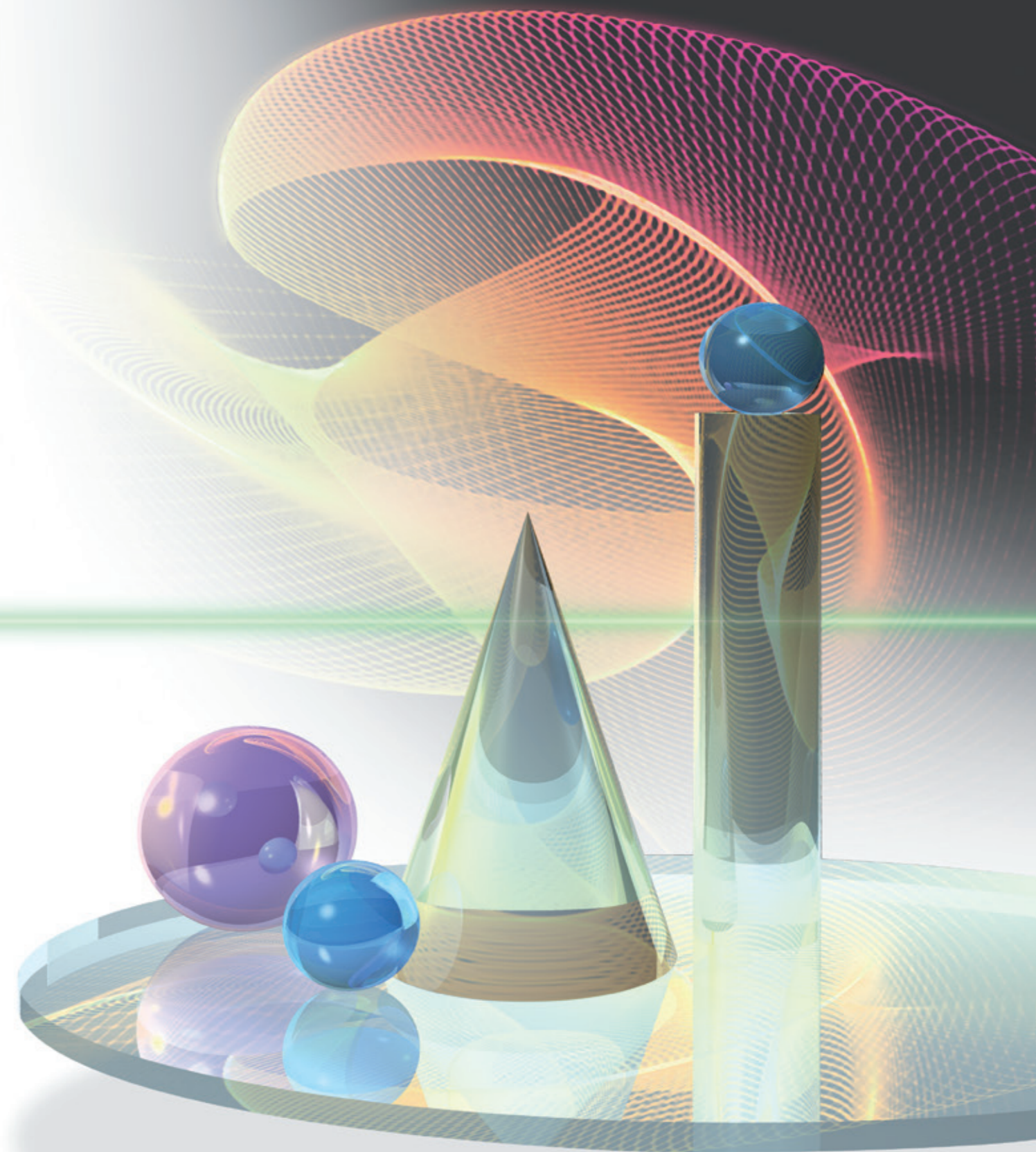Meiritz Seiki has been specializing in the manufacturing of vibration isolation systems since 1971. Ever since, we have designed vibration isolators to protect precision instruments from the harmful effects of vibration. Our focus has always been to develop new technologies that meet future market demands—a philosophy that has helped us continue maintain our expertise and leadership in this field. This is particularly true for optical experiment applications, where Meiritz was the first company in Japan to utilize the aluminum honeycomb design. To help us better meet the diversifying needs of our customers, we have recently complemented our product line with a series of steel honeycomb tables. Included in this catalog is the entire line of our products intended for use in basic optical experiments. We feel confident that the catalog will enable you to find the right equipment for your needs.

### CONTENTS

除振 Vibration Isolation	3
空気ばね式除振装置 Air Spring Vibration Isolators	4
ハニカム定盤 Honeycomb Surface Table	5
AYA Series	7
AY/AYN Series	11
RHS/LHS Series	15
HS/HSB Series	21
AET Series	24
K6 Series	27
G Series	28
関連機器	29

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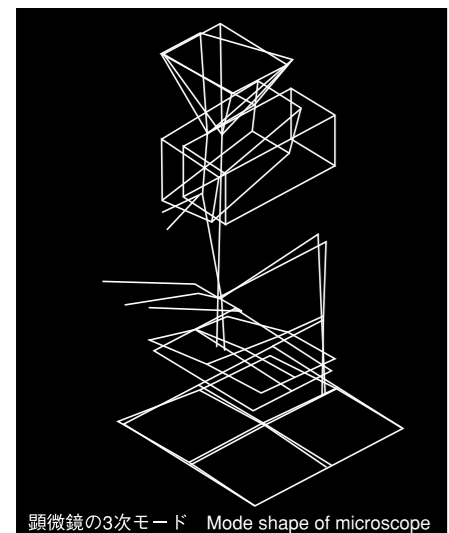
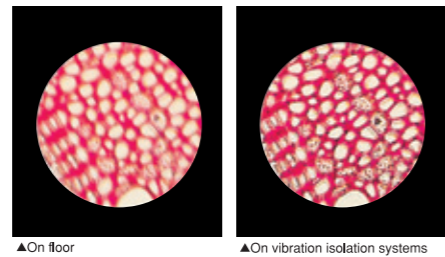
The most effective means of dealing with vibrations for precision instruments is to interrupt the vibrations immediately prior to the precision instrument itself.

What is Harmful Vibration in the case of Precision Instruments?

Precision instruments are composed of numerous components each having their own natural resonance frequency. If vibration at the resonance frequency of a particular component is transmitted to that component, that component begins to vibrate excessively causing a disturbance in the positional relationship of all of the other components. For example, when an analysis of the vibration of a metal microscope was performed, it was confirmed that the microscope had a particular mode in which the lens barrel of the microscope curved slightly at a frequency of 130Hz. Therefore, the resonance frequency of the lens barrel is 130Hz, and vibration at this resonance frequency can be considered to be harmful vibration which must be isolated from the microscope.

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
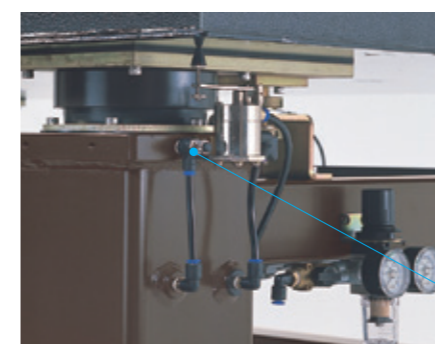
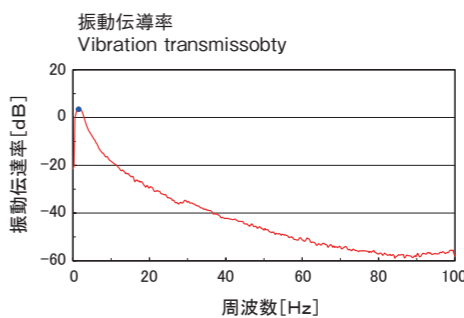
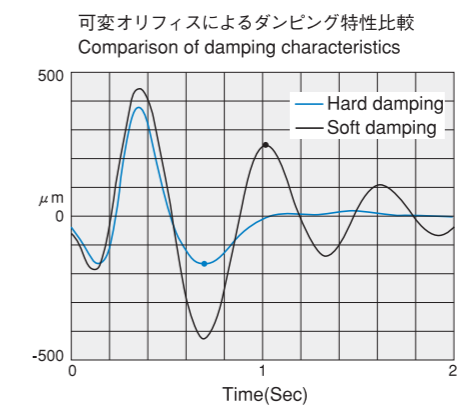
Air Spring System Featuring Leading-Edge Technology

In determining the performance of an air spring system, factors such as the shape and material of the air springs, the volume of the air spring chambers and the surge tank, the diameter and length of the orifice, and the type of automatic leveler are all mutually inter-related. Therefore, a high level of technical capability and a vast amount of experience are necessary in realizing superior performance. All of the isolation system models house the finest in air spring systems that can only be made possible through the expertise of Meiritz. allows it to respond with a high degree of sensitivity to even extremely light objects, minimizing vibration spiraling as much as possible.

Patented Automatic Leveler

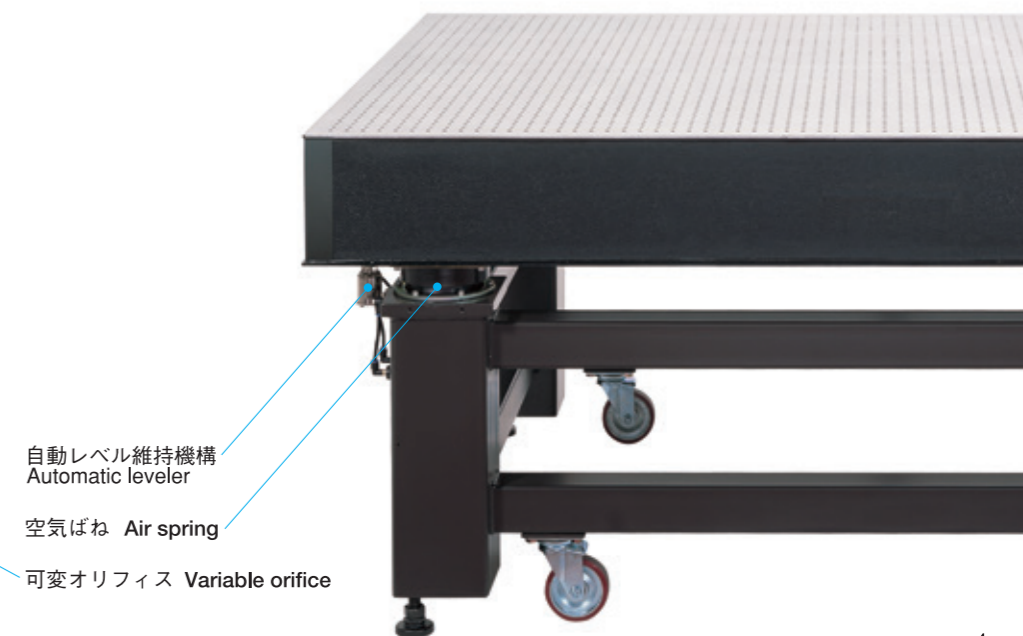This series of vibration isolators are equipped with our patented automatic leveling mechanism which constantly keeps the surface table level even when removing and placing instruments as well as when changing their position. The unmatched performance of this automatic leveler allows it to respond with a high degree of sensitivity to even extremely light objects, minimizing vibration spiraling as much as possible.

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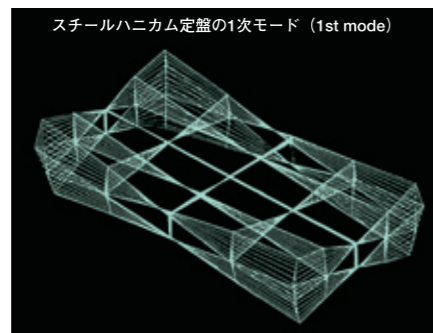
### Highly-Advanced Steel Honeycomb Table

Following overall improvement of the damping materials and the steel honeycomb cores, our steel honeycomb table has evolved into a highperformance optical table that offers extreme stiffness and damping. Reflecting our environmental concerns, a bonding agent having no organic solvents is used between the honeycomb cores and the top and bottom steel plates.

### Machine Polishing Offers Maximum Table Surface Flatness

Machine polishing of a honeycomb table surface, which has been regarded as difficult to perform, has become possible thanks to our unique technology. As an optional feature, tables with maximum flatness can be manufactured to order. (please contact us for more information).

### We Meet Your Diversified Application Requirements

The tabletop length, width and thickness can be modified upon request. Work surface options and configurations, such as through holes, cut-outs, notches, and custom tapped holes, are also available. Two or more tables can also be joined together to create longer or wider tables, or T-shaped tables, L-shaped tables, etc. We have manufactured a variety of complex tables in different shapes, on request.

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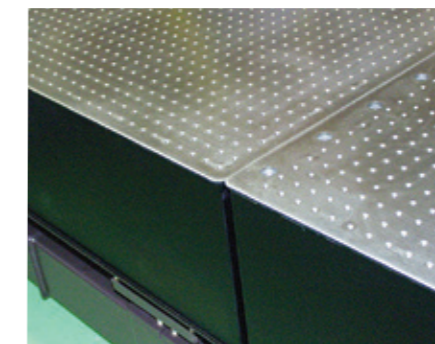
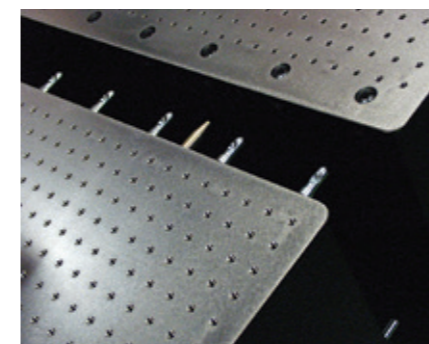
### Type K4

This high-performance model is the result of many years of experience and innovation. The top plate is made of SUS430 stainless steel, enabling the use of a magnetic stand. Type K4 comes standard with M6 tapped holes for mounting optical equipment, either on 50mm grid or 25mm grid (to be specified by the user).

### Type K5

Type K5 is based on Type K4. Each M6 tapped hole, either on 50 mm grid or 25mm grid(to be specified by the user), is sealed with a cap attached on the back side, to prevent inflow of water or chemicals, and the circulation of air by convection into the honeycomb table. The air spring vibration isolator of Type K5 (AYA and AY series) comes standard with an automatic leveler with exhaust air drain line, for use in clean rooms.

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The AYA Series is Meiritz's top-of-the-line series which features improved isolation performance and various useful functions.

The AYA Series is the crystallization of advanced technology developed entirely by Meiritz. In particular, through improvement of horizontal direction vibration isolation performance, the AYA Series is able to respond to the optical experimentation needs including interferometers and holography equipment which are sensitive to minute vibrations in all directions.

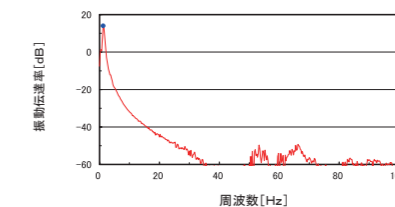
### Parallel Swing Mechanism

The AYA Series is equipped with Meiritz's original parallel swing mechanism which can be likened to a swing which can swing forward and backward as well as to the right and left, and moreover, can do so simultaneously. The result is the realization of the natural frequency in the horizontal direction being reduced to 0.5 Hz, less than 1/3 that of our previous models.


### New Mechanisms for Emphasis on Ease of Use

The AYA Series is equipped with a newtype of damper which utilizes the new material αGEL Damping performance, which is able to promptly absorb any vibrations which cause the surface table to fluctuate, has been improved even more without any loss of vibration isolation performance thereby leading to increased ease of use. αGEL is the registered trademark of Taika Co.,Ltd

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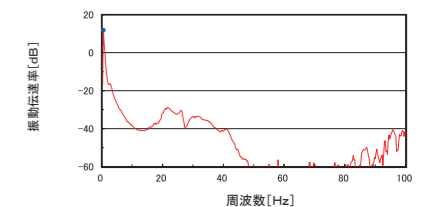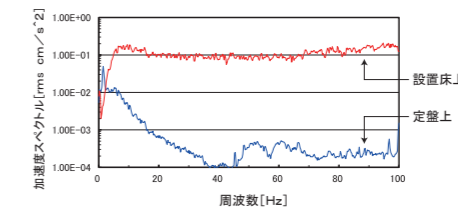
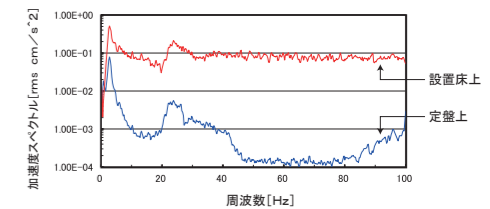
#### AYA-Type K4

- M6 tapped holes on 25 mm grid are manufactured as standard. (Can also manufacture 50 mm grid)
- Through hole and special tapping work available. (Optional)
- Can join honeycomb tables with no deterioration of stiffness. (Optional)
- Can use a magnet stand.

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#### AYA-Type K5

- Caps are attached to the back of M6 tapped holes to seal the holes. Equipped with the concentrated exhaust type of automatic leveler sensor as standard.
- M6 tapped holes on 25 mm grid are manufactured as standard. (Can also manufacture 50 mm grid)
- Can join honeycomb tables with no deterioration of stiffness. (Optional)
- Can use a magnet stand.

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This popular series gives top priority to cost performance. Meiritz has made this series of vibration isolators in order to provide an inexpensive vibration isolator that can easily be used in the field. Such development concept has been realized through setting specifications in consideration of standardization and mass production, as well as thorough rationalization of processing and assembly processes.

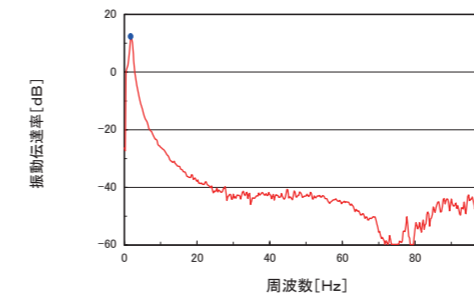
### Unmatched Reliability Supported by Leading-Edge Technology

Meiritz makes a careful check of various performance conditions taking full advantage of their vast amount of accumulated data relating to floor vibrations and vibration of equipment placed on vibration isolator, as well as vibration analysis technology. This has allowed us to succeed in standardizing the specifications that are truly required by the numerous vibration isolation needs of today. As a result, we have realized the providing of vibration isolators that feature high reliability plus a reasonable price.


### Manual Leveling Types also Available

There are certain users who require large vibration isolators for use in rooms without air sources. In response to this demand, Meiritz has made available a series of vibration isolators that allow leveling of the surface table by supplying air with an attached pump.

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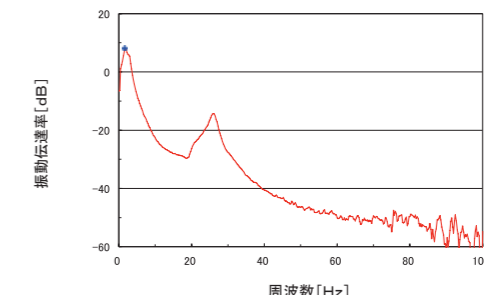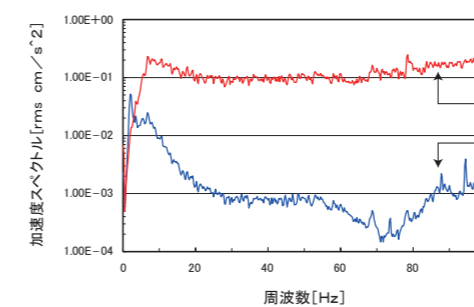
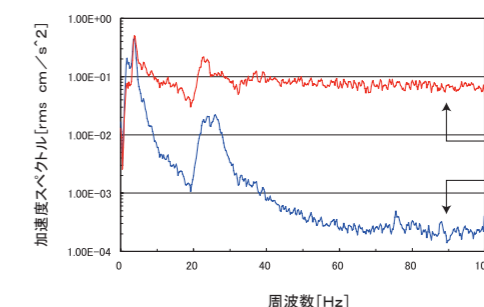
#### AY/AYN-Type K5

- Caps are attached to the back of M6 tapped holes to seal the holes. Equipped with the concentrated exhaust type of automatic leveler sensor as standard.
- M6 tapped holes on 25 mm grid are manufactured as standard.  
(Can also manufacture 50 mm grid)
- Can join honeycomb tables with no deterioration of stiffness. (Optional)
- Can use a magnet stand.

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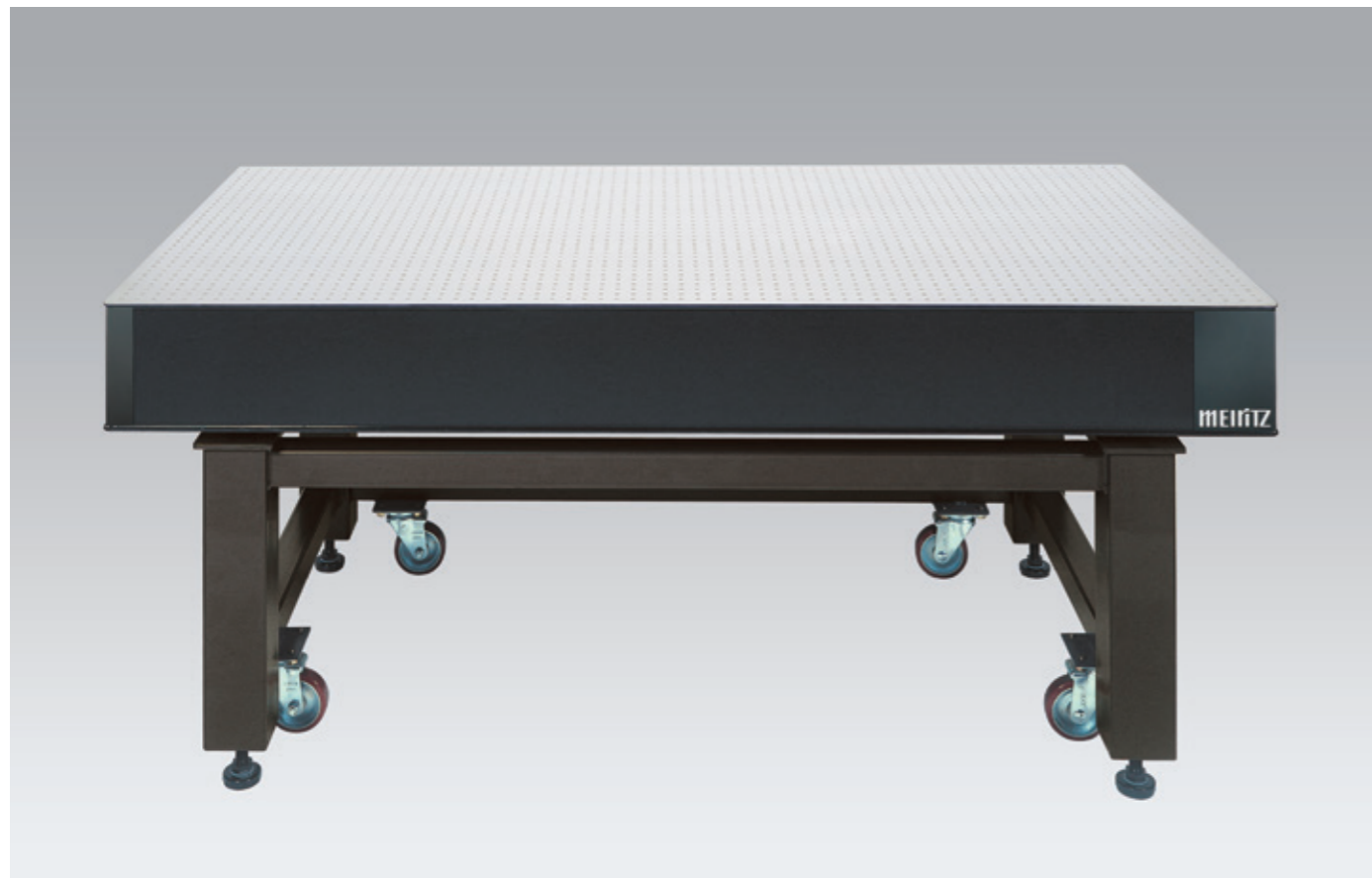
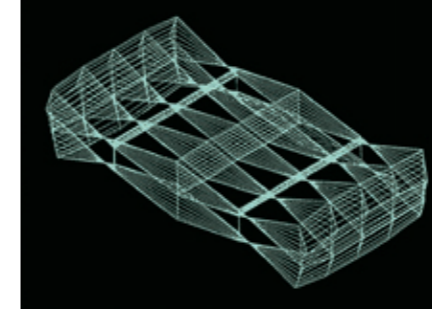
These surface tables specially designed for optical experimentation applications have inherited all of the characteristics of the surface tables used in our steel honeycomb surface table air spring vibration isolators, equipped with all of the features found in these long-running, best-sellers.

### RHS Series-Optical Surface Tables with Stand

The stand provided with these optical surface tables utilizes vibration analysis technology cultivated in the design of vibration isolators. It features an optimal design that brings out the features of optical surface tables. Since the stand is equipped with foot-adjusted and casters, the installation site can be changed easily.

The RHS Series uses a rubber cushion to suppress vibrations. Rigid tables without rubber cushions, called the LHS Series, are also available.

### Layout markings and Tap Processing Available (Optional)

Inscription of marking lines useful in determining the layout of the equipment to be placed on the surface table, as well as tap processing for anchoring of equipment are available on an optional basis (M5, M6)

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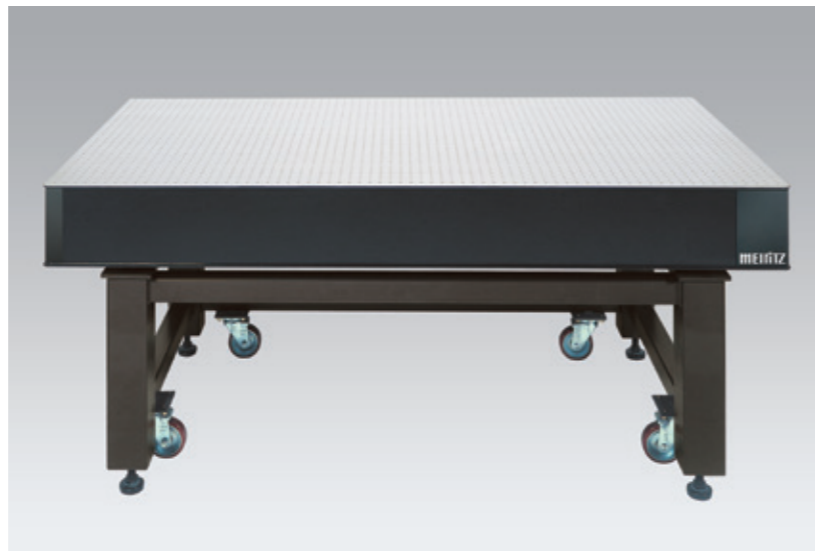
#### LHS-Type K4 (Rigid type without rubber cushion)

- M6 tapped holes on 25 mm grid are manufactured as standard.  
(Can also manufacture 50 mm grid)
- Through hole and special tapping work available. (Optional)
- Can join honeycomb tables with no deterioration of stiffness. (Optional)
- Can use a magnet stand.

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#### LHS-Type K5 (Rigid type without rubber cushion)

- Caps are attached to the back of M6 tapped holes to seal the holes.
- M6 tapped holes on 25 mm grid are manufactured as standard.  
(Can also manufacture 50 mm grid)
- Can join honeycomb tables with no deterioration of stiffness. (Optional)
- Can use a magnet stand.

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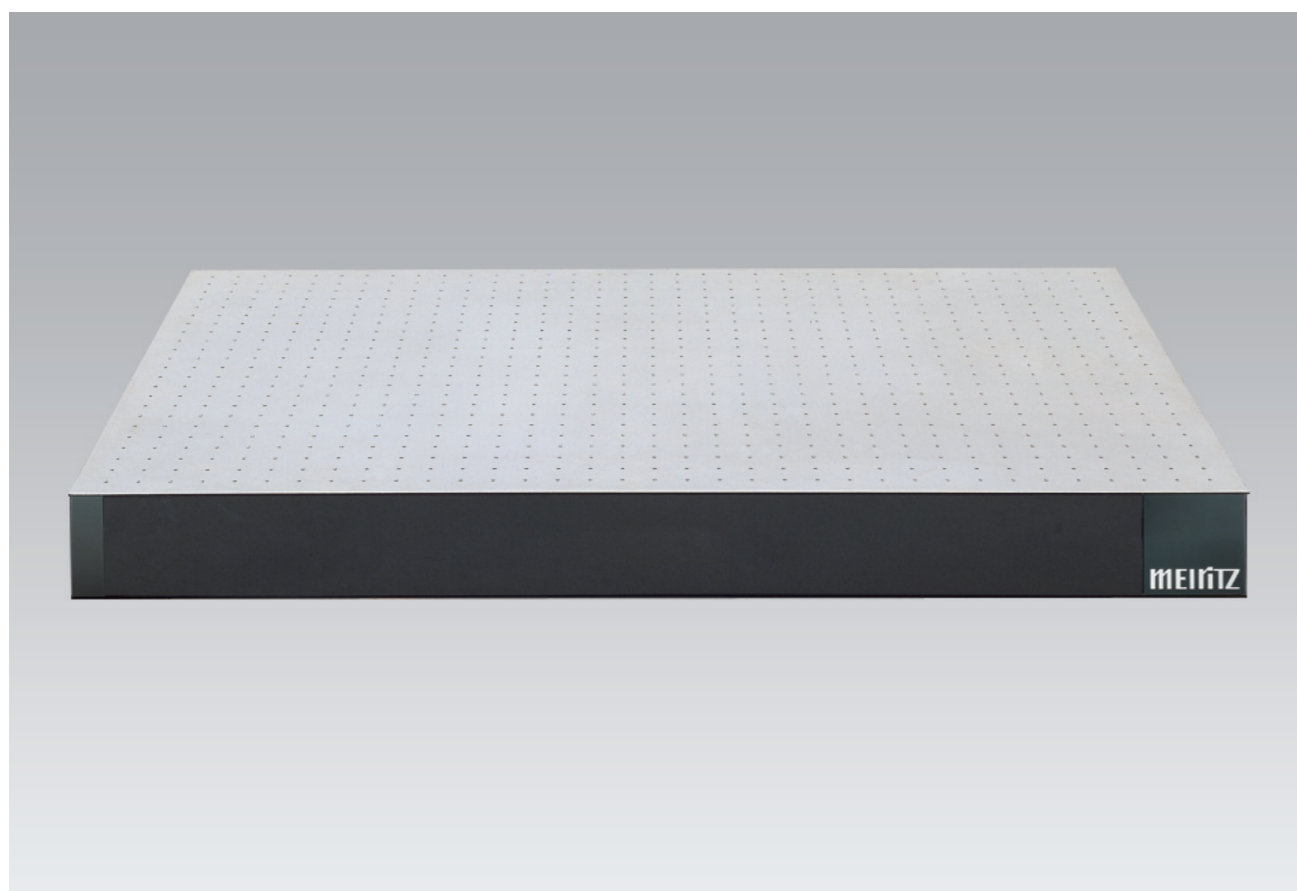
HS-Type K4

- M6 tapped holes on 25 mm grid are manufactured as standard (Can also manufacture 50 mm grid)
- Through hole and special tapping work available. (Optional)
- Can join honeycomb tables with no deterioration of stiffness. (Optional)
- Can use a magnet stand.
- 4 rubber plates are attached.

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RHS-Type K5

- Caps are attached to the back of M6 tapped holes to seal the holes.
- M6 tapped holes on 25 mm grid are manufactured as standard. (Can also manufacture 50 mm grid)
- Can join honeycomb tables with no deterioration of stiffness. (Optional)
- Can use a magnet stand.
- 4 rubber plates are attached.

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
HSB-Type K4

- M6 tapped holes on 25 mm grid are manufactured as standard. (Can also manufacture 50 mm grid)
- Through hole and special tapping work available. (Optional)
- Can use a magnet stand.
- 4 rubber plates are attached.
- M6 tapped holes can be sealed by caps. (Optional)

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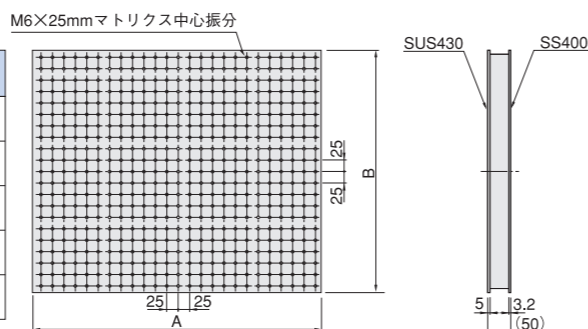The welding structure stand has been added to the thin steel honeycomb table, a time-proven model from the HSB Series, to achieve low cost and high stiffness.

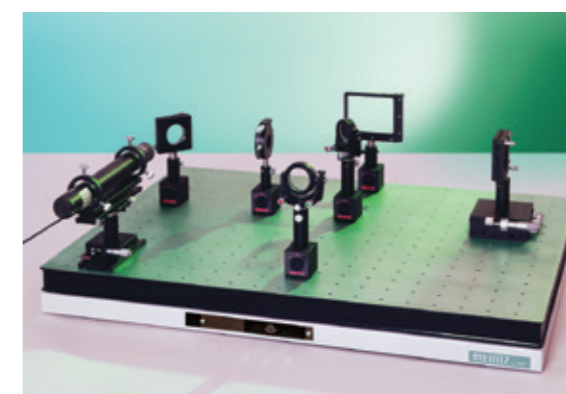
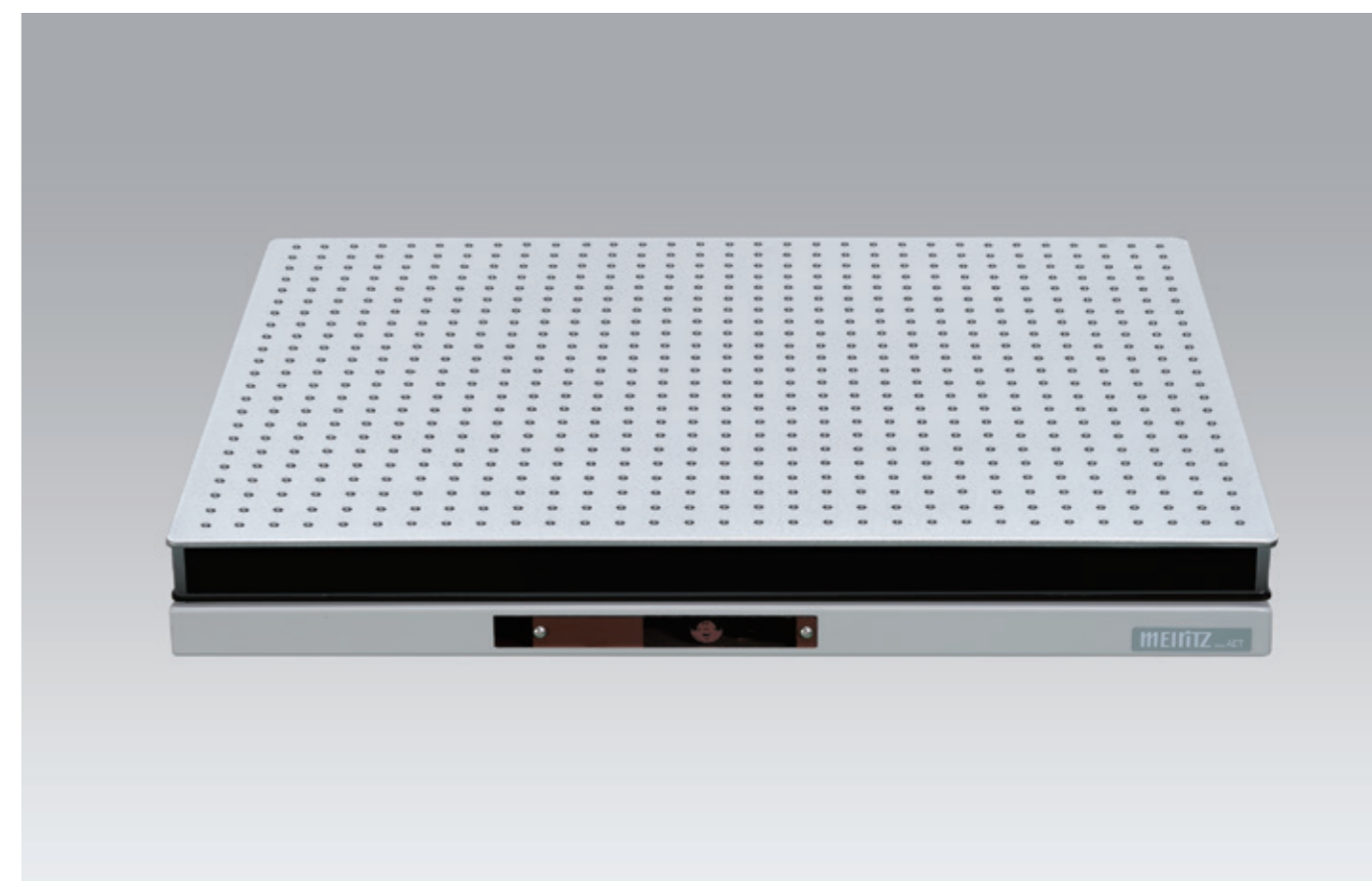
※Through holes with special specifications can be manufactured on request. Please contact us for available dimensions and specifications.

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Business Fields

- For Microscopy
- For Optical & Laser
- For Measurement

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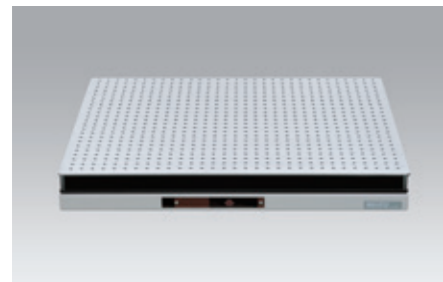
The AET Series is a series of desktop vibration isolators that bring together an entire high-performance vibration isolation system. At Meiritz Seiki, many of our previous desktop vibration isolators have been combined into a series. As a result, the AET Series, which combines these vibration isolators with steel honeycomb surfaces tables, consists of desktop air suspension vibration isolators are optimal for optical experiments using a small surface area.

## Four Types Available

### ● S Type with Automatic Leveler

The S type is equipped with an automatic leveler. This keeps the surface table level even when the load distribution on the spring system is changed. The automatic leveler is constantly connected to an air source of 0.35-0.70Mpa and the air pressure can be reduced with the pressure control valve provided.

### ● Manual Leveling N Type

The N type is provided with a hand pump used for supplying air to the system in order to maintain the surface table level. We recommend this type for use in locations not equipped with an air source.

### ● Clamp-Equipped NC and SC Type

The SC and NC types are equipped with clamps that allow you to fix the surface table in position both quickly and easily.

### ● H Type with high load specifications

Specifications of a load weight of up to 200 kg have been added to the lineup to deal with heavier devices. (Size 0405 can load up to 100 kg)

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The other critical element for an optical table, besides high stiffness, is high damping performance. In other words, it is important to damp the vibrations as soon as they are received by the table. Meiritz Seiki offers a new optical table that offers even higher damping performance than its tried and tested steel honeycomb table.

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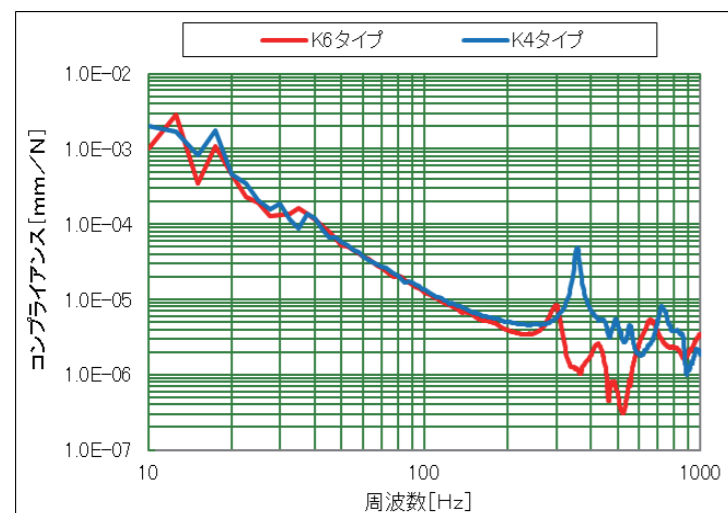
Outline and features

- high stiffness, is high damping performance. In other words, it is important to damp the vibrations as soon as they are received by the table. Meiritz Seiki offers a new optical table that offers even higher damping performance than its tried and tested steel honeycomb table.

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
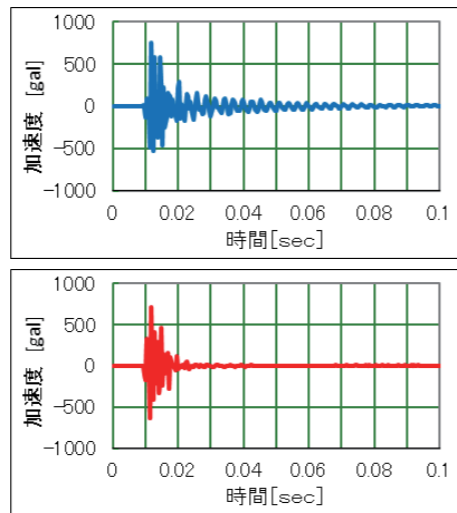
Flexible combination with stands

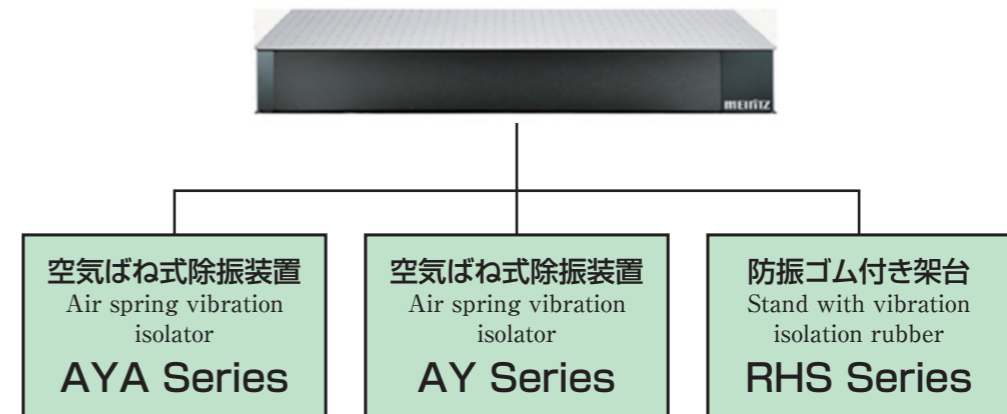The Steel Honeycomb Table K6 Type enables flexible combination with a stand equipped with an air spring vibration isolation mechanism from the AYA and AY Series or with an RHS stand with vibration isolation rubbers.  
(※Tables of the same size as the K4 Type or of special sizes can be manufactured on request.)

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The granite table offers outstanding features such as high surface accuracy and little secular distortion and, moreover, it is barely influenced by changes in environmental conditions such as temperature and humidity. It is suitable for precision machinery and precision measurement that places particular emphasis on the mechanical and physical characteristics of the table.

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Main features of granite table

- High surface accuracy is obtained.
- The thermal conductivity is low, and there is little impact from surrounding temperature changes.
- The absorption rate is low and the table is not susceptible to humidity.
- Even if the table is scratched by mistake, the surface will not become burred and the surface flatness will be maintained.
- Granite is rarely worn out because it is twice or two and a half times as hard as cast iron.
- The elastic coefficient is high, offering excellent deflection stiffness.

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**MEIRITZ SEIKI CO.,LTD.**